



ETHANOL GAS STANDARD CYLINDER REPORT
 NORTH DAKOTA OFFICE OF ATTORNEY GENERAL
 CRIME LABORATORY DIVISION-TOXICOLOGY SECTION/BREATH ALCOHOL PROGRAM
 SFN 59282 (08/2018)

Chemical Test Operator Name (Print) <i>Patrick Thomas</i>		
Location <i>Fargo Police Department</i>	Intoxilyzer® Serial Number <i>80-004942</i>	
Gas Lot Number <i>19825080A2</i>	Gas Cylinder Number <i>33</i>	Gas Expiration Date <i>09/05/2027</i>

Check When Done:

- 1. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- AEI 2. Perform an ACA Test (Level 1, Function C).
 02Mar26 Write in the ACA Test Results (Reported to 3 Digits, Ex. 0.081).
 - A. 0.080 AC
 - B. 0.079 AC
 - C. 0.080 AC
- 3. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.
- 4. File Previous Intoxilyzer® Record (SFN50496, Form 120-G) at the Intoxilyzer® Location at the Agency.
- 5. Send the Following to the Crime Laboratory:
 - A. Ethanol Gas Standard Cylinder Report (SFN59282, Form 105-G).
 - B. ACA Test Record.

Chemical Test Operator Signature <i>Patrick Thomas</i>	Date <i>02/25/2026</i>
Reviewed By (Crime Laboratory Use Only) Anna Ingemansen <i>Anna Ingemansen</i>	Date 02March2026

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-004942
Location = FAPD 8164.14.00 09/16
02/25/2026 17:07

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	17:08
02 Std. Gas	0.080	17:09
03 Room Air	0.000	17:09
04 Std. Gas	0.079	17:09
05 Room Air	0.000	17:10
06 Std. Gas	0.080	17:10
07 Room Air	0.000	17:11

Lot No = 19825080A2
Cyl No = 33
Exp Date = 09/05/2027
County = 09 Oper No. = 130056


Operator Signature
PATRICK THOMAS

Remarks: *ACA all within acceptable ranges*

Form 106-I8000