



INTOXILYZER® 8000 INSTALLATION AND REPAIR CHECKOUT
 NORTH DAKOTA OFFICE OF ATTORNEY GENERAL
 CRIME LABORATORY DIVISION-TOXICOLOGY SECTION/BREATH ALCOHOL PROGRAM
 SFN 59281 (06/2018)

Serial Number 80-005937	Instrument Location Belcourt PD
Reason for Install/Repair <input checked="" type="checkbox"/> Install After Receiving From Crime Laboratory <input type="checkbox"/> Install After Location Change <input type="checkbox"/> Other (Specify) _____	

Check When Done:

- 1. Surge Protector Installed/Property Grounded.
- 2. Telephone Line Connected to Intoxilyzer® 8000.
- 3. Breath Tube Heated.
- 4. Enter Preliminary Data (i.e. Date, Time, DST (Y), and Location; Level 2, Function E).
- 5. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- 6. Run Tests:
 - A. Print Test (Level1, Function P).
 - B. ACA Test (Level 1, Function C).
 - C. Radio Frequency Interference (RFI) Test (CMS Mode or Level 1, Function B or C; Key Radio During Test).
- 7. Repair and/or Maintenance Performed (if any): _____
- 8. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.
- 9. File Previous Intoxilyzer® Record (SFN504096, Form 120-G) at the Intoxilyzer® Location at the Agency.
- 10. Send the Following to the Crime Laboratory: Completed Intoxilyzer® 8000 Installation and Repair Checkout (SFN59281, Form 104-G), Print Test, ACA Test, and RFI Test.

Field Inspector Signature <i>Jerrid Marin</i> - Jerrid Marin	Date 4/8/26
-----------------------------------------------------------------	-----------------------

Crime Laboratory Use Only

This installation has been reviewed and the instrument is approved to be used for the analysis of breath to determine alcohol concentration from the date the Field Inspector performed the installation. This record on file at the Office of Attorney General, Crime Laboratory Division, in the County of Burleigh, North Dakota, is certified to be a true and correct copy of the documents received.

Reviewed/Certified By Anna Ingemansen <i>Anna Ingemansen</i>	Certified Date 01May2026
------------------------------------------------------------------------	------------------------------------

The initial print test, ACA Test and RFI Test had a typo in the Standard Lot #. This was corrected and a new print test is attached to reflect the update. This does not affect the installation of this instrument. AEI 01May2026

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005937
Location = BELC 8164.14.00 09/16
04/08/2026 11:05

***** Printer Test *****

abcdefghijklmnopqrstuvwxyz1234567890-=
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#%^&*()_+?

abcdefghijklmnopqrstuvwxyz1234567890-=
ABCDEFGHIJKLMNPOQRSTUVWXYZ!@#%^&*()_+?

Current Instrument Setup


Data Entry Mode: Enabled
Start Test Sequence: DABACABA
Display Prelim Rslt? Yes
Display Third Digit? Yes
Inhib Printer(Y/N)? No
Display Volume? No
Disable On Memfull? Yes
of Print Copies? 1
Select Std (D/W/I)? Dry
Standard Value? 0.080
Standard Lot #? 08425008A1
Standard Cyl #? 20
Standard Expiration? 05/05/2027
Oper No? 080929

Flow Cal. Date: 11/17/2017
Slope 606
Intercept -534563

IR Calibration Date: 11/17/2017
 3um 9um

0th Coef(*100): -30534 -23772
1st Coef(*100): 294336 142357
2nd Coef(*100): -906 850
H2O adj(mg/l*10k): 205 250

***** Printer Test End *****

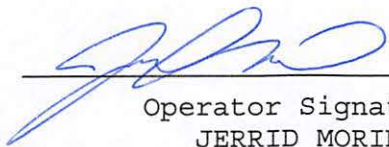

Operator Signature
JERRID MORIN

Remarks:


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005937
Location = BELC 8164.14.00 09/16
04/08/2026 11:05

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	11:06
02 Std. Gas	0.079	11:07
03 Room Air	0.000	11:07
04 Std. Gas	0.079	11:07
05 Room Air	0.000	11:08
06 Std. Gas	0.080	11:08
07 Room Air	0.000	11:09

Lot No = 08425008A1
Cyl No = 20
Exp Date = 05/05/2027
County = 63 Oper No. = 080929


Operator Signature
JERRID MORIN

Remarks:

Form 106-I8000


Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-005937
Location = BELC 8164.14.00 09/16
04/08/2026 11:09

Test	AC	Time
01 Room Air	0.000	11:11
02 *Subject Test	RFI*	11:11
03 Room Air	0.000	11:12

*Invalid Test
Inhibited - RFI

Sub Name = TEST, TEST NA
Sub DOB = 01/01/1999
Sub Sex = Male Weight = 250
Test = DUI Cit = NA
Dr. Lic. = ND/POI000000
Lot No = 08425008A1
Cyl No = 20
Expiration Date = 05/05/2027
County = 63 Oper No. = 080929

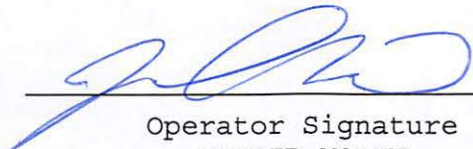

Operator Signature
JERRID MORIN

Remarks:

Form 106-I8000

