



ETHANOL GAS STANDARD CYLINDER REPORT

NORTH DAKOTA OFFICE OF ATTORNEY GENERAL

CRIME LABORATORY DIVISION-TOXICOLOGY SECTION/BREATH ALCOHOL PROGRAM

SFN 59282 (08/2018)

Chemical Test Operator Name (Print) <u>Howard, Austin L</u>		
Location <u>Burke county Sheriff's Shop</u>	Intoxilyzer® Serial Number <u>80-006496</u>	
Gas Lot Number <u>264437</u>	Gas Cylinder Number <u>118</u>	Gas Expiration Date <u>1/29/28</u>

Check When Done:

- ☒ 1. Scan/Enter Gas Cylinder Information (Level 1, Function S).
- ☒ 2. Perform an ACA Test (Level 1, Function C).
Write in the ACA Test Results (Reported to 3 Digits, Ex. 0.081).
 - A. 0.082 AC
 - B. 0.079 AC
 - C. 0.079 AC
- ☒ 3. Complete the Top Portion of the Intoxilyzer® Record (SFN50496, Form 120-G) and Place it by the Intoxilyzer® for Use.
- ☒ 4. File Previous Intoxilyzer® Record (SFN50496, Form 120-G) at the Intoxilyzer® Location at the Agency.
- ☒ 5. Send the Following to the Crime Laboratory:
 - A. Ethanol Gas Standard Cylinder Report (SFN59282, Form 105-G).
 - B. ACA Test Record.

Chemical Test Operator Signature <u>Austin Howard</u>	Date <u>6/5/26</u>
Reviewed By (Crime Laboratory Use Only) <u>Anna Ingemansen</u> <u>Anna Ingemansen</u>	Date <u>17 Jun 2026</u>

Intoxilyzer Test Record and Checklist
NDOAG Crime Lab. Div., Bismarck, ND 58501

CMI, Inc. Intoxilyzer Alcohol Analyzer
North Dakota Model 8000 SN 80-006496
Location = BURK 8164.14.00 09/16
06/05/2026 17:07

DRY CAL CHECK

Test	AC	Time
01 Room Air	0.000	17:08
02 Std. Gas	0.080	17:08
03 Room Air	0.000	17:09
04 Std. Gas	0.079	17:09
05 Room Air	0.000	17:10
06 Std. Gas	0.079	17:10
07 Room Air	0.000	17:11

Lot No = 264437
Cyl No = 118
Exp Date = 01/29/2028
County = 07 Oper No. = 132545



Operator Signature
AUSTIN HOWARD

Remarks:

Form 106-I8000